OLM-056 Stabförmige LED-Strahlungsquelle Rod-shaped LED radiant source

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The OLM-056 Module is a rod-shaped LED radiation source with an integrated battery. The module contains six UV-LEDs. Each LED has a glass lens that focus the light at an angle of about 20°. For an effective cooling an aluminum heat sink is used underneath the LED-plate. Pressing the power button on the top will turn on the LED Module and allow you to irradiate the surface area you aim at for a desired time. This easy-to-use LED Module is the best possible solution for various applications in research, development and industrial applications, for example:

- Phototherapy and other medical applications
- Disinfection of surfaces
- Florescence excitation
- Photo and bio-chemistry

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The system is designed for a fast implementing and usage in laboratory or for medical issues.

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This System contains the following parts

- LED module OLM-056
- Power supply for the battery
- Connecting cable

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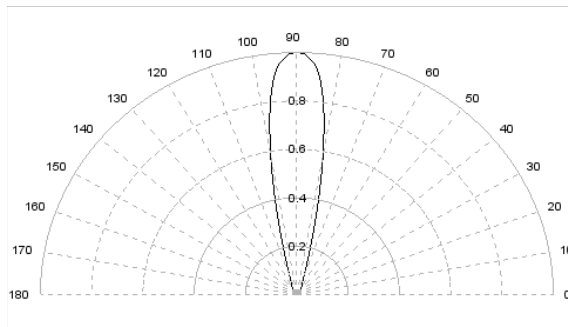
- Size 262 x 40 x 41 mm
- Material: aluminum, anodized
- Light emitting area: 75 x 27 mm
- LED-number: 6
- Heat sink material: aluminum
- Weight: about 500 g
- Window: quartz glass

Type	Peak wavelength ¹⁾ (nm)	Typical power density ²⁾ (mW/cm ²)	Max current per LED Chip (A)	Forward Voltage (V)
OLM-056D265	260 – 270	15	0.5	39
OLM-056D310	305-315	20	0.5	36
OLM-056D365	360-370	200	0.5	21
OLM-056D400	395-405	300	0,5	20


¹⁾ Other wavelength in the range 265-1050 nm can be supplied upon request

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
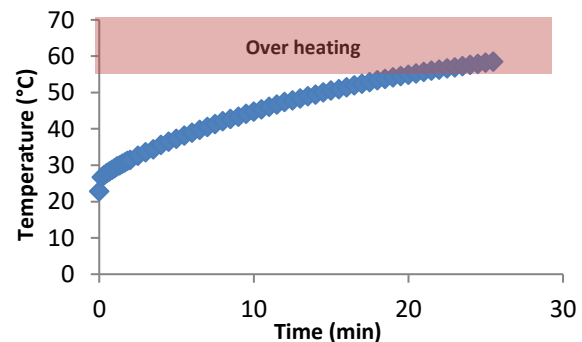
Temperature development @ 365 nm

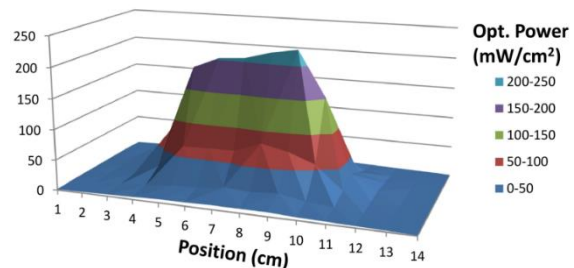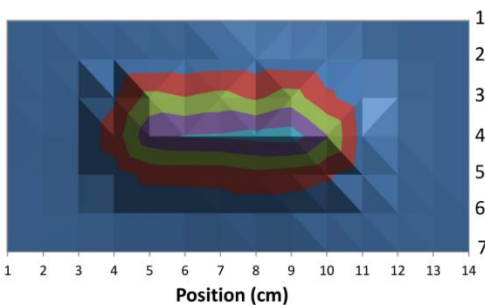The LED Module contains an internal temperature sensor. In combination with the control electronics the module is switched off at 55°C till the temperature drop below 45°C and therefore protected against over temperature

Optical power distribution in a distance of 3 cm to the glas plate @ 365 nm

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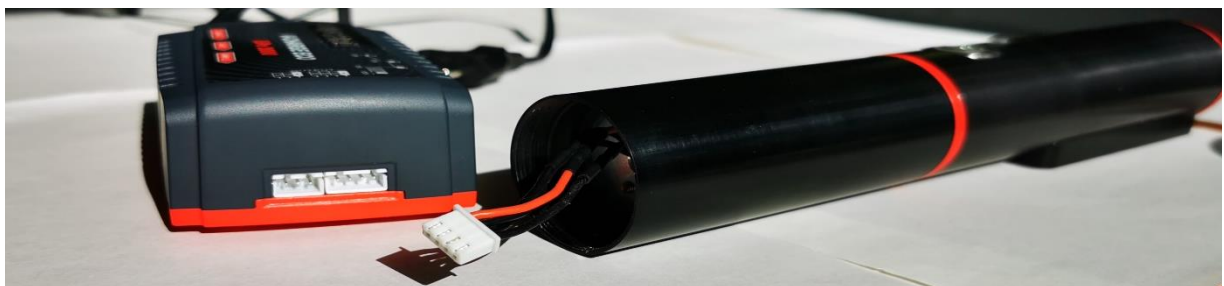
- 1) Quartz glass (75 x 27 mm)
- 2) 6x LED with lens (365 nm)
- 3) Aluminum housings
- 4) Battery with cable for charging
- 5) Control LED-display when ON/OFF Button pressed:
 - Green (permanent): System functioning well
 - Yellow (blinking): battery almost empty
 - Yellow (permanent): battery empty
 - Red (permanent): overheating
- 6) Power button (ON / OFF) with ring illumination (white)

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- Voltcraft V-Charge Eco LiPo 2000
- Dimensions: 110 x 70 x 40 mm
- weight: 1610g
- Operating voltage: 100-240V / AC
- Charging current (max.): 2A
- Charging capacity: 25 W
- Short-circuit protection

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- Capacity: 1000 mAh
- Power rating: 25C
- Weight: 104 g
- Connecting system: BEC
- Balancer-Connecting system: XH

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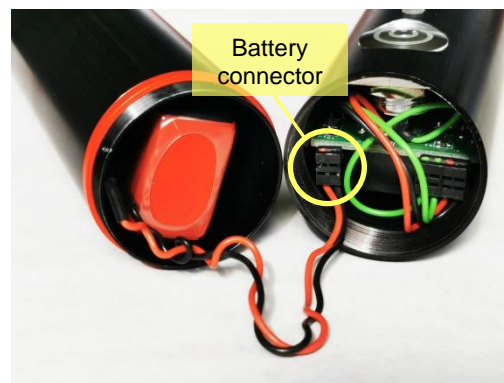
Bringing into service

- Read carefully the operating manual and the safety instructions
- The press button (6) turns module on / off. The ring illumination indicates the operation of the module
- At over temperature the module will be switched off and the control LED-display (5) will turn red.
- The Control LED-display (5) will blink yellow to indicate that the battery should be recharged. In case the LED-Display is permanently yellow, the battery is empty and has to be charged immediately to prevent deep discharge.

Shutting down

After operation if the LED Module is not going to be used in the next 24 hours, disconnect the battery from the control electronic plate to avoid deep discharge. Fail to do so will make the battery in a few days out of order without a possibility for repair.

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- These Power LEDs emit during operation intense UV optical radiation, which may be harmful to eyes, even for brief periods.
- Do not look directly into the LED during operation
- Be sure that you and all persons in the vicinity wear suitable eye protection, face mask and skin protection (e.g. gloves) if the radiation of the LED is not covered by other measures
- LED are not a toy, keep out of reach of children
- The user has the responsibility to inform, train and instruct customers and co-workers

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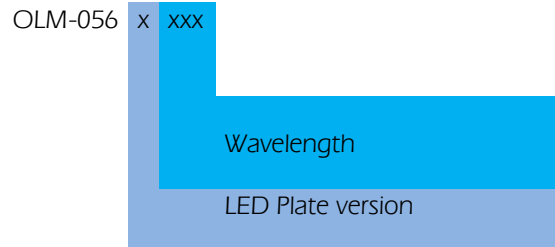
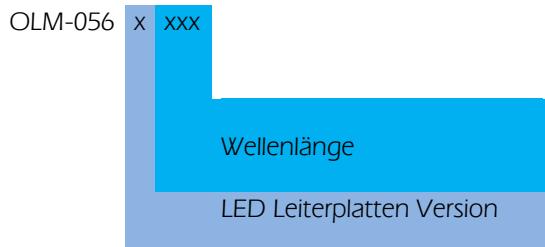
www.osa-opto.com Made in Germany

CE

CAUTION
**Ultraviolet LED
 Radiation**
 Do NOT expose to
 unprotected eyes or skin